



Technical Bulletin

FMCG Industry Solutions Pty Ltd ABN 22 135 446 007 Factory 11A, 1 - 3 Endeavour Rd, Caringbah NSW 2229 www.fmcgis.com.au sales@fmcgis.com.au 1300 628 104 or 02 9540 2288

Delvotest® Fast BT

Sensitivity of Delvotest® Fast BT

Delvotest® Fast BT is test for qualitative detection of antibiotic residues in raw milk, namely \(\text{B-lactams} \) and tetracyclines. Delvotest® Fast BT is based on immunochromatographic technology to qualitatively detect antibiotic residues in milk. The list below is not exhaustive and valid for Delvotest® Fast BT, the sensitivity of the test is given only for the most commonly used antibiotics. The level of detection is based on 100% positive.

Raw Cow Milk

Antibiotic class	Antibiotic	S/TL (US)	MRL (EU)	Level of detection Cow
		ppb		
Penicillins	Penicillin G	5	4	2-3
	Ampicillin	10	4	3
	Amoxicillin	10	4	3
	Oxacillin	-	30	10
	Dicloxacillin	-	30	10
	Nafcillin	-	30	15
	Cloxacillin		30	5
Cephalosporins	Ceftiofur	50	100	4
	Cefquinome	-	20	10
	Cefazoline	-	50	25
	Cephapirine	20	60	10
	Cefacetrile	-	125	50
	Cefoperazone	-	50	10
	Cefalexine	-	100	1000
	Cefalonium	-	20	4
Tetracyclines	Tetracycline	300	100	60-80
	Oxytetracycline	-	100	80
	Chlortetracycline	-	100	>200
	Doxycycline	-	0	>1000

Although diligent care has been used to ensure that the information provided herein is accurate, nothing contained herein can be construed to imply any representation or warranty for which we assume legal responsibility, including without limitation any warranties as to the accuracy, currency or completeness of this information or of non-infringement of third party intellectual property rights. The content of this document is subject to change without further notice. This document is non-controlled and will not be automatically replaced when changed. Please contact us for the latest version of this document or for further information. Since the user's product formulations, specific use applications and conditions of use are beyond our control, we make no warranty or representation regarding the results which may be obtained by the user. It shall be the responsibility of the user to determine the suitability of our products for the user's specific purposes and the legal status for the user's intended use of our products.

DSM Food Specialties B.V. | A. Fleminglaan 1 | 2613 AX Delft | The Netherlands | info.food@dsm.com | Trade Register Number 27235314





Raw Goat and Sheep milk

Antibiotic class	Antibiotic	MRL (EU)	Level of detection Goat	Level of detection Sheep	
		ppb			
Penicilins	Penicillin G	4	1.5	2	
	Ampicillin	4	3	3	
	Amoxicillin	4	3	3	
	Oxacillin	30	10	nd	
	Cloxacillin	30	5	5	
Cephalosporins	Cefazoline	50	25	15	
	Cephapirine	60	10	10	
Tetracyclines	Tetracycline	100	40	80	
	Oxytetracycline	100	40	100	

Although diligent care has been used to ensure that the information provided herein is accurate, nothing contained herein can be construed to imply any representation or warranty for which we assume legal responsibility, including without limitation any warranties as to the accuracy, currency or completeness of this information or of non-infringement of third party intellectual property rights. The content of this document is subject to change without further notice. This document is non-controlled and will not be automatically replaced when changed. Please contact us for the latest version of this document or for further information. Since the user's product formulations, specific use applications and conditions of use are beyond our control, we make no warranty or representation regarding the results which may be obtained by the user. It shall be the responsibility of the user to determine the suitability of our products for the user's specific purposes and the legal status for the user's intended use of our products.

DSM Food Specialties B.V. | A. Fleminglaan 1 | 2613 AX Delft | The Netherlands | info.food@dsm.com | Trade Register Number 27235314