

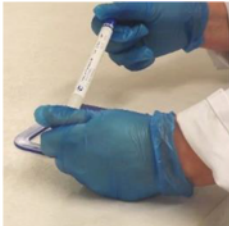

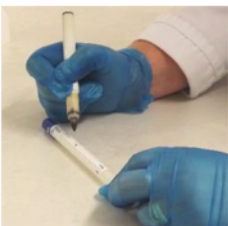
GUIDANCE FOR USE

SwabSURE Ecoli/Coliform – SS-L04



FMCG Industry Solutions Pty Ltd
 ABN 22 135 446 007
 Factory 11A, 1 – 3 Endeavour Rd, Caringbah NSW 2229
 www.fmcgisl.com.au
 sales@fmcgisl.com.au
 1300 628 104 or 02 9540 2288

All samples should be kept cool until testing is performed. Please visit the website for Precautions/Limitations.

<p>STEP 1</p> <p>Identify the site to be sampled. It is recommended to use a template for accurate results where possible (TS/15-T).</p> 	<p>STEP 2</p> <p>Hold the swab with both hands, pull down and twist the swab cap away from the tube; breaking the security seal in the process.</p> 
<p>STEP 3</p> <p>Remove swab from tube using the cap as a holder. Do not touch the swab tip or shaft and make sure that it does not touch any surface before sampling. Do not discard the tube.</p> 	<p>STEP 4</p> <p>Sample designated area using a template where possible. If the area permits, swab in 3 planes (horizontal, vertical and diagonal) whilst rotating swab between forefinger and thumb.</p> 
<p>STEP 5</p> <p>When sampling is completed insert swab back into the tube and push cap on tightly.</p> 	<p>STEP 6 (DELAYED TEST)</p> <p>Write sample details on tube label. Store at 4-8°C (or cool & dry as possible) and send to the laboratory within 24 hours to initiate the chromogenic test.</p> 
<p>STEP 7 (IMMEDIATE TEST)</p> <p>Alternatively (where permitted), immediately insert swab into desired test media tube until the breakpoint is level with the top of the vial. Pull down on the swab shaft whilst holding the tube straight to activate the breakpoint. Replace Cap & agitate lightly.</p> 	<p>STEP 8 (CHROMOGENIC TEST)</p> <p>Place tubes into a rack & incubate them vertically at 36±1°C.</p> 